

HiPette

Fully Autoclavable Mechanical Pipette

The new generation of DLAB fully autoclavable manual pipette is lightweight, precise, stable and durable. It adopts new piston structure, metal screw structure, fluorine rubber compound sealing ring and other technologies to reduce the pipetting resistance. At the same time, it can withstand high temperature and high pressure sterilization, resist strong chemical corrosion, and have a long service life. In addition, volume lock helps achieve more precise pipetting operations.







Complete sterilization

Strong materials ensure HiPette can fully autoclavable, UV sterilizable and resistent strong chemical corrosion.



Forward pipetting

Suitable for low-viscosity liquids, press the plunger to the first stop when aspirating, and press the button to the second stop when dispensing.

Reverse pipetting

It is suitable for high-viscosity liquids, easy to foam or micro samples, with higher pipetting accuracy. Press the plunger to the second stop when aspirating, and press the plunger to the first stop when dispensing.

Range selection

Choose the appropriate range according to the pipetting volume. 30%—100% of the maximum range of the pipette is the best use range.

Specifications

Single-channel Adjustable Volume Pipettes (9 models, covering 0.1μl~10ml.)						
Volume Range	Increment	Test Volume	Accuracy Error		Precision Error	
			%	μL	%	μL
0.1~2.5μL	0.002μL	2.5µL	1.4%	0.04	0.7%	0.02
		1.25µL	2.5%	0.03	1.5%	0.02
		0.25μL	12.0%	0.03	6.0%	0.02
0.5~10μL	0.01μL	10μL	1.0%	0.10	0.4%	0.04
		5μL	1.5%	0.08	0.8%	0.04
		1μL	2.5%	0.03	1.8%	0.02
2~20μL	0.02μL	20μL	1.0%	0.20	0.3%	0.06
		10μL	1.2%	0.12	0.6%	0.06
		2μL	5.0%	0.10	1.5%	0.03
10~100μL	0.1μL	100μL	0.8%	0.80	0.2%	0.20
		50μL	1.0%	0.50	0.3%	0.15
		10μL	3.0%	0.30	1.0%	0.10
20~200μL	0.2μL	200μL	0.6%	1.20	0.2%	0.40
		100μL	1.0%	1.00	0.3%	0.30
		20μL	2.5%	0.50	0.7%	0.14
30~300μL	2μL	1000μL	0.6%	1.80	0.2%	0.60
		500μL	1.0%	1.50	0.3%	0.45
		100μL	2.5%	0.75	0.7%	0.21
100~1000μL	1μL	1000μL	0.6%	6.00	0.2%	2.00
		500μL	1%	5.00	0.2%	1.00
		100μL	3%	3.00	0.6%	0.60
1000~5000μL	5μL	5000μL	0.6%	30.00	0.2%	10.00
		2500μL	0.6%	15.00	0.3%	7.50
		1000μL	1.2%	12.00	0.3%	3.00
2000~10000μL	10μL	10000µL	0.6%	60.00	0.2%	20.00
		5000μL	0.8%	40.00	0.2%	10.00
		2000μL	1.5%	30.00	0.3%	6.00

^{*} For accurate and reliable pipette calibration HiPette conforms to ISO 8655:2002 standards.